

General remarks

The Cytometry Facility location in room WAD-K-106 is a BL-2 working area. Food and drinks are not allowed inside the rooms! All waste containers must be decontaminated and neutralized before disposing of any liquids.

Solid Waste

- Solid BSL-1 and BSL-2 waste is collected in the red waste container
- Do not fill the red waste containers with non-biohazardous waste
- Non-biohazardous waste can be disposed of in the standard trash bins.
Any trash placed in a standard trash bin must be safe for the cleaning personnel.

Aria cell sorter (liquid waste)

- **Decontaminated waste containers must be treated before disposal!**

Last user of the day at Aria cell sorter

- Disconnect full waste container from the instrument and attach the empty waste container to the instrument
- Add the appropriate volume of Bleach/Javel, (located under the sink) to the tank
- Put the magnetic “bleach added” label on the tank.
- Store the tank close to the sink in WAD-K-106 for overnight incubation.

LSR Fortessa (liquid waste)

- **Decontaminated waste containers must be treated by the facility staff before disposal!**

User changing the Fortessa Waste upon alarm goes off on the wet cart

- Disconnect full waste container from the instrument and attach the empty waste container to the instrument.
- Add 100 ml of Bleach/Javel, (located under the sink) to the tank.
- Put the magnetic “bleach added” label on the tank.
- Store the tank close to the sink in WAD-K-106 for overnight incubation.

Cytek Aurora (liquid waste)

- **Decontaminated waste containers must be treated by the facility staff before disposal!**

User changing the waste tank when prompted by the software

- Disconnect full waste container from the instrument and attach an empty box to the instrument
- Add **200 ml** of Na-hypochlorite/Bleach/Javel (blue container located under the sink).
- Write a note on the tank that bleach has been added.
- Store the tank close to the sink in WAD-K-106 for overnight incubation.

Amounts of Na-hypochlorite/Bleach/Javel to add:

Tank	Fill level	Amount of bleach to add
10 l plastic tanks (BD instruments)	¼ full	25 ml
	½ full	50 ml
	full	100 ml
20 l carton box (Aurora)	full	200 ml